



Vector® Fusion-Aid™ - HA Kit
Catalog No. MB-0734

This Fusion-Aid™ kit is designed to isolate fusion proteins containing HA-tag. This kit contains an affinity-purified antibody to HA-tag coupled through a stable hydrophilic spacer arm to crosslinked 4% agarose beads. Spin columns and collection tubes are included for convenience. For optimal results it is recommended that the protocol on the reverse side of this specification sheet be followed.

Product **AGAROSE ANTI-HA**
AFFINITY PURIFIED

Amount 0.5 ml of settled gel (1:1 v/v slurry)

Lot No. ZA0520

Produced in goat

Suspension Solution 10 mM phosphate, 0.15 M NaCl, pH 7.5, 0.08% sodium azide.

Antibody Concentration 2.0 mg/ml of settled gel

Binding Capacity ≥0.4 mg HA tagged β-galactosidase (120kD)/ml

Storage Conditions Refrigerate - DO NOT FREEZE

Other Remarks

*Actual binding capacity may vary depending on the size of the fusion protein.

Related Products

	<u>Cat. No.</u>	<u>Unit Size</u>
Biotinylated Anti-HA, affinity purified, made in goat	BA-0704	0.25 mg
VECTASTAIN® ABC-AmP™ Reagent (Standard Kit)	AK-6000	1 kit
DuoLUX™ Chemiluminescent/Fluorescent Substrate for Alkaline Phosphatase	SK-6605	100 ml
BCIP/NBT Substrate Kit	SK-5400	1 kit
10x Casein Solution	SP-5020	250 ml
Biotinylated Protein Molecular Weight Markers	SP-1400	50 blots